


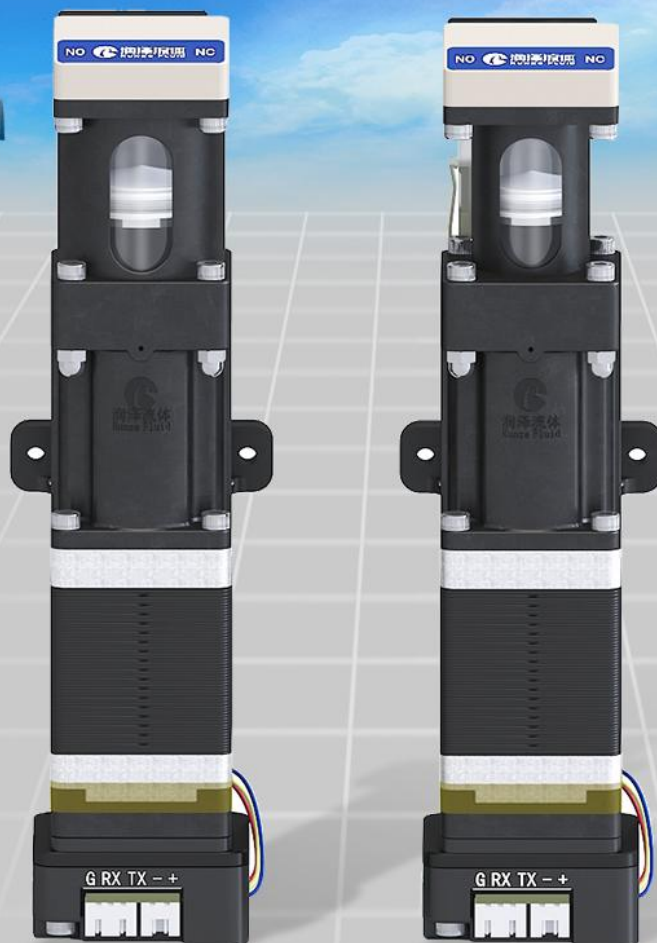

V 1.2

SY-09S Vertical Syringe Pump Manual

南京润泽流体控制设备有限公司
NANJING RUNZE FLUID CONTROL EQUIPMENT CO.,LTD

Note: This manual applies to RUNZE protocol. If you are not sure of the current pump using the protocol, please check P20[C query and jump protocol] first

Table of Contents

Chapter 1 product introduction	2
1.1 Company Overview	3
1.2 SY-09S Features at-a-glance	3
1.3 Product Safety Precautions	3
Chapter 2 product introduction	4
2.1 Product Features	5
2.2 Naming Rules	5
2.3 Basic Parameters	5
2.3.1 Product Function	5
2.3.2 Technical Parameters	6
Chapter 3 hardware setting	7
3.1 Component Installation	8
3.1.1 Structure Diagram	8
3.2 Manual for 35 Stepper Motor	8
3.3 Port Definition	10
Chapter 4 software communication	11
4.1 Brief Introduction	11
4.2 Installation & Debugging	11
4.3 Communication Protocols	11
4.3.1 Command format	11
Chapter 5 common problems & solutions	17
Chapter 6 quick command	18
A Quick Commands	18
B Hexadecimal Quick Reference Table	19
C Switching Protocol	20
Chapter 7 version description	22
Chapter 8 technical service	23

Chapter 1 Product Introduction

1.1 Company Overview

Nanjing Runze Fluid Control Equipment Co., Ltd., established in 2014, is a national high-tech enterprise focusing on R&D and production of fluid accessories for numerous analytical instruments. We engineer, manufacture and market differentiated standard products such as syringe pumps, multiport valves, peristaltic pumps, gastight syringes, plastic fittings, etc. We persevere in providing our customers with best quality and service in the fields of environmental monitoring, biopharmaceuticals, medical equipment, industrial automation and laboratory instruments, etc.

In past years, we have accumulated rich technical and practical experience that bring us honors of ISO9001, National High-tech Enterprise, Jiangsu Province Private Science and Technology Enterprise, 5A Bank Credit Assessment, 44 technical patents and software copyright including 2 invention patents, 27 utility model patents, 13 design patents, 2 software copyright.

1.2 SY-09S Features at-a-glance

Congratulations on your purchase of the SY-09S Syringe Pump from RUNZE Fluid.

SY-09S Syringe Pump is a fully programmable, small compact size, high-precision liquid handling micro industrial pump module with stable performance & long service life, developed by RUNZE Company. Controlled by a host controlling system (external computer, microprocessor, PLC, etc.), the clockwise or counterclockwise circular motion of the stepper motor is converted into linear motion through the trapezoidal screw rod, which makes the syringe pump piston move up and down linearly to achieve aspirating and dispensing functions.

Configuration : 3ml, 8ml

Component : Borosilicate glass syringe, trapezoidal screw, optocoupler, stepper motor, drive.

Usage : SY-09S syringe pump is widely used in liquid transferring system with high-precision and high-stability sampling requirements, such as laboratory instrument, medical analysis equipment, chromatographic analyzer, automatic biochemical analyzer, blood analyzer, trace element analyzer, electrolytic analyzer, food & beverages detection and analysis system, water quality on-line analyzer, petroleum detection equipment and biopharmaceutical extraction devices.


1.3 Product Safety Precautions

1. The device needs to be reset with 0x4F for the next command after power on; if any abnormality occurs during the use, please also reset with 0x4F command.

2. Applicable power supply: DC 24V $\pm$ 10%, 3A, when using linear power supply, the voltage and current

must be adjusted to the corresponding parameter value

3. When debugging this product by computer, please adopt RUNZE debugging software serialcomm
4. In the debugging process, the liquid must be debugged to avoid dry wear of the piston, affecting the service life of the injector
5. When using this product, please connect the ground wire to reduce the interference caused by environmental factors.
6. Please use the original serial cable of this product to connect with the power supply.
7. The two communication modes of this product (RS232 and RS485 bus) are in non-isolated mode.
8. Do not disassemble the product parts, tamper-proof label torn without warranty.
9. When operating the software, please refer to the software operating instructions and communication protocols, and do not make up data input without permission.
10. Please dispose of the instrument in accordance with the regulations for disposal of instrument and equipment waste. Users should not throw it away at will.
11. When connecting multiple devices with RS485 bus protocol, please refer to the connection method in Figure 1-1 below, but the resistance value should be decided according to the number of devices hooked up by the user.

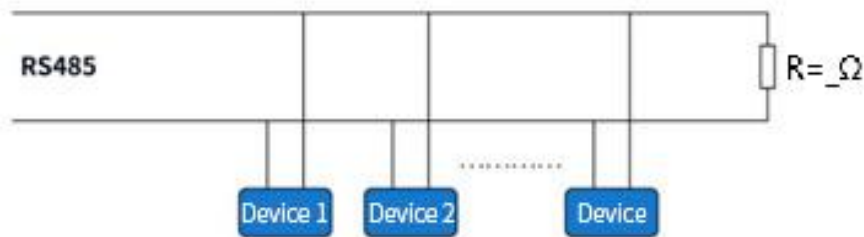

Chapter 2 Product Introduction

2.1 Product Features


Small size, space-saving installation, high precision, stable performance, long service life of micro industrial syringe pump

The contact material is made of high borosilicate glass and PTFE, which is corrosion resistant and high temperature resistant and suitable for many special media

Widely used in environmental protection equipment, medical analysis equipment, non-standard high-precision sampling equipment and other analyzers

2.2 Naming Rules

Parameters

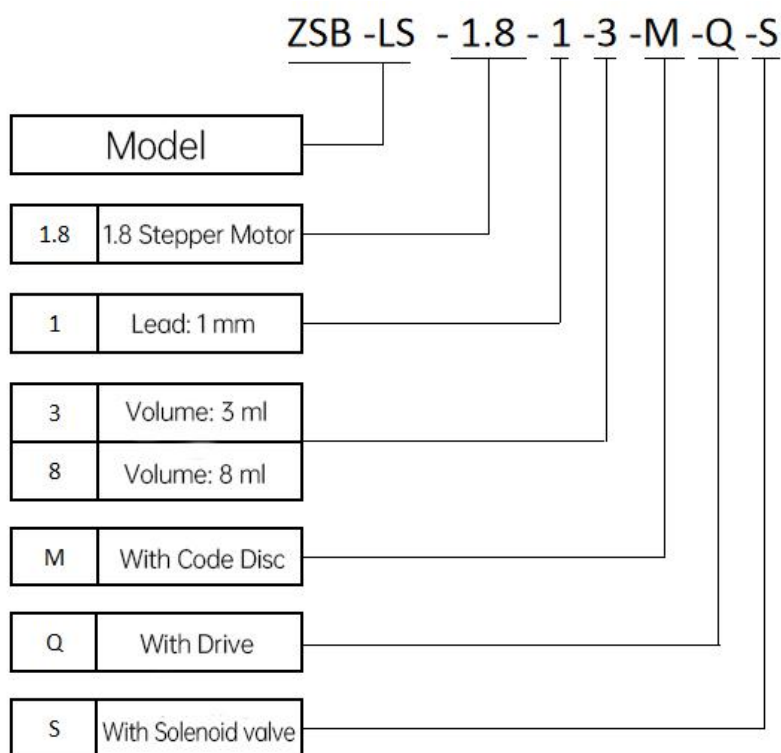

For example: the 3ml syringe pump, with 1.8 degree stepper motor, single hole, female thread with driver is named ZSB-LS-1.8-1-3-M-Q

2.3 Basic Parameters

2.3.1 Product Function

Product Function	Description
Set address	Set serial port address.
Set baud rate	Different baud rates for RS232, RS485

Set CAN destination address	If there are multiple devices, the CAN address can be set as the priority of different devices.
Set speed	3ml : from 1rpm to 600rpm 8ml : from 1rpm to 300rpm (There are difference for gas, liquid and models.)
Set subdivision	Subdivision can be set from subdivision 2 to 32.
Reset internal data	Restore factory settings.
Query parameter	The device address, speed, subdivision and baud rate can be queried.
Query version	Query the current firmware version.
Motor control direction	The motor can be controlled to rotate clockwise and counterclockwise.
Reset	Return the piston of the syringe pump to the home position.
Force stop	Stop the current operation of the syringe pump motor.
Query motor status	Check the current motor status.

2.3.2 Technical Parameters


Product Function	Description	
Accuracy	≤1% (rated stroke)	
Precision	0.3%-0.7% (rated stroke)	
Service life	3 million times no leakage (media: water; 1 rated stroke = one time)	
Volume	3ml	8ml
Rated stroke (Control steps)	18mm(3600 步)	19.2mm(3840 步)
Maximum speed	600rpm	300rpm
Linear speed	0.017~10mm/s	0.017~5mm/s
Running time (Per rated stroke)	1.8~1080s	3.84~1152s
Resolution	0.005mm/0.833μl	0.005mm/2.083μl
Syringe ID	14.55mm	23.03mm
Actuator	Trapezoidal screw (Lead 1mm)	
Wetted material	Borosilicate glass, PTFE, PPS, PEEK, EPDM, FKM	
Maximum pressure	0.075~0.2	
Channel	Single channel	
Connection	1/4-28UNF	
Communication interface	RS232/RS485	
Baud rate	RS232/RS485 : 9600bps/19200bps/38400bps/57600bps/115200bps	
Address & Parameter	Via communication	

setting		
Power supply	DC24V/3A	
Rated power	15W	
Operating temperature	5 ~ 55C°	
Operating humidity	≤80% (relatively humidity, non-condensing)	
Dimension (L*W*H)	51*45.8*177.4	51*45.8*176.21
Weight	0.56kg	0.62kg


Chapter 3 Hardware Setting

3.1 Component Installation

3.1.1 Structure Diagram

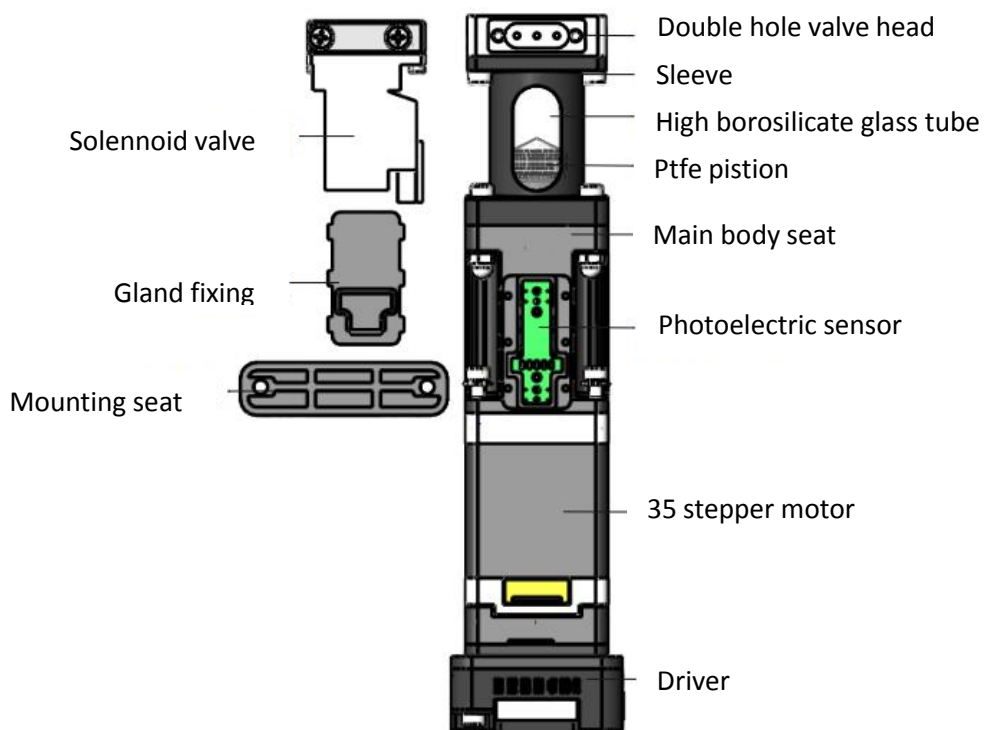

3.1.2 Dimension without Driver (Unit: mm)

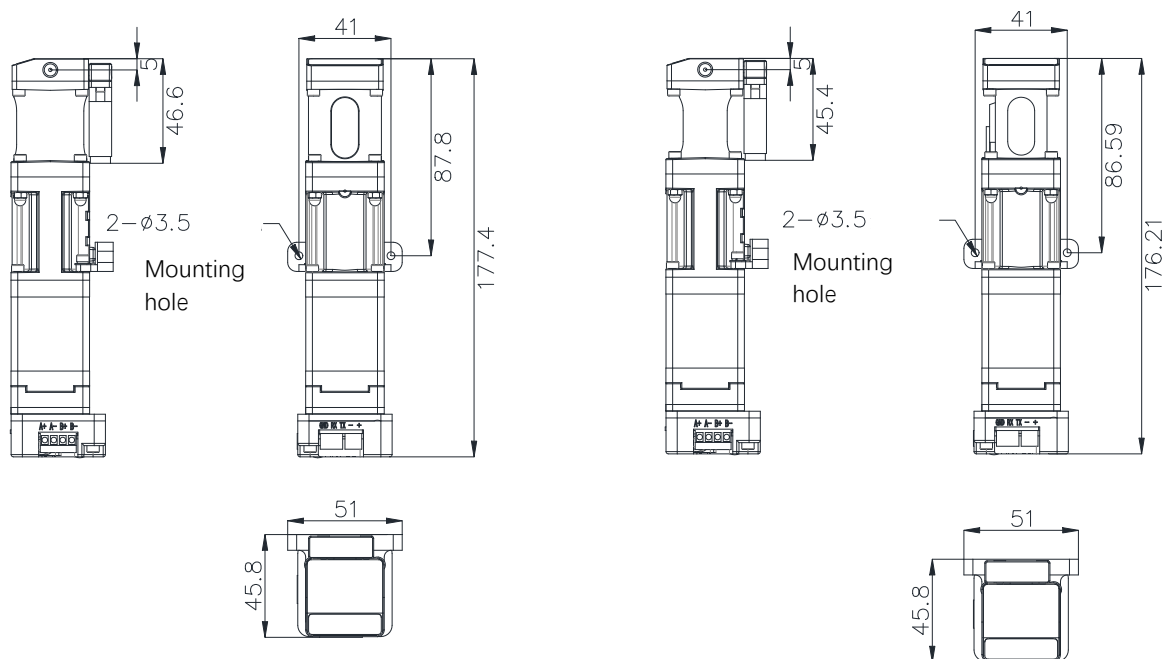

3.2 Manual for 35 Stepper Motor

L/W STEP	赤 RED	青 BLU	黄 YEL	橙 ORG
1	+	+	-	-
2	-	+	+	-
3	-	-	+	+
4	+	-	-	+

FIG. 2


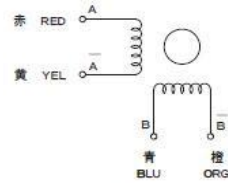


FIG. 1

出 版
I S S U E D
'20.11.03
STEPPING MOTOR
MinebeaMitsumi Inc.

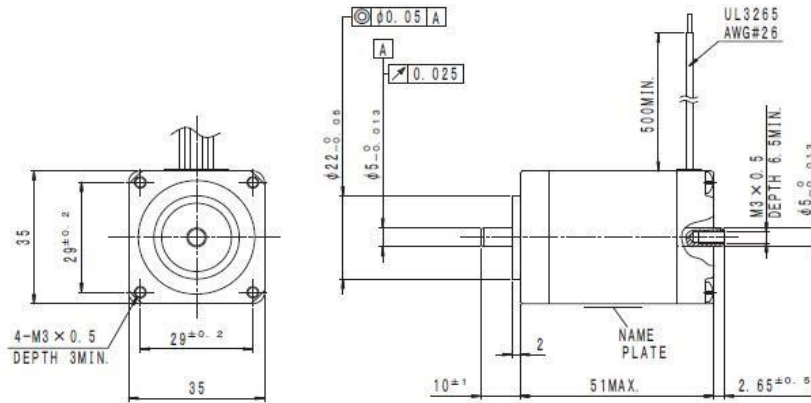

ミネベアミツミ株式会社 MinebeaMitsumi Inc.	UNIT	材質 MATERIAL	作成日 DATE
APPROVED	検査 CHECKED	検査者名 NAME	OCT. / 30 / 20
IGRS	SATO	検査場所 PLACE	品名 DESCRIPTION
PHOL	TEL	検査設備 EQUIPMENT	HYBRID STEP MOTOR
		検査結果 RESULT	図番 DRAWING NO.
			14PM-F447CNWA5607

6. 信頼性 Reliability


下記条件で試験後、2項の電気的特性、3項の機械的特性を満足するものとする。

The product to be examined in the following condition and satisfy 2.Electrical characteristics and 3.Mechanical characteristics.

No.	試験項目 Item	試験条件 Test condition	参考規格 Reference standard
1	低温放置 Cold (耐寒性)	温度 Temperature : $-40^{\circ}\text{C} \pm 3^{\circ}\text{C}$ 試験時間 Test time : 96 h	JIS C 60068-2-1
2	高温放置 Dry heat (耐熱性)	温度 Temperature : $85^{\circ}\text{C} \pm 2^{\circ}\text{C}$ 試験時間 Test time : 96 h	JIS C 60068-2-2
3	高温高湿放置 Damp heat (耐湿性)	温度 Temperature : $60^{\circ}\text{C} \pm 2^{\circ}\text{C}$ ($40^{\circ}\text{C} \pm 2^{\circ}\text{C}$) 湿度 Humidity : 93 % +2 / -3 % 試験時間 Test time : 96 h	JIS C 60068-2-3
4	冷熱衝撃 Thermal shock (温度変化)	温度 Temperature : 低温側 Cold $-40^{\circ}\text{C} \pm 3^{\circ}\text{C}$ 1 h ↑ ↓ (温度移行時間: 5 min 以内) (Temperature migration time: 5 min MAX.) 高温側 Dry heat $85^{\circ}\text{C} \pm 2^{\circ}\text{C}$ 1 h サイクル数 Number of cycle : 25 cycle	JIS C 60068-2-14
5	振動 Vibration	振動数範囲 Frequency range : 10 Hz ~ 55 Hz ~ 10 Hz 振 幅 Amplitude : 1.5 mm 掃 引 時 間 Time coefficient: 1 min 加 振 方 向 Direction of excitation: X, Y, Z 時 間 Time : 各方向 Each direction 2 h	JIS C 60068-2-6
6	衝撃 Shock	加 速 度 Acceleration : 981 m/s ² { 100 G } 正弦半波 Half-sine 作 用 時 間 Interaction time: 6 ms 落 下 方 向 Drop direction: $\pm X, \pm Y, \pm Z$ 落 下 回 数 Number of drop: 各3回 Each 3 times	JIS C 60068-2-27

3.3 Port Definition

Diagram of the driver control board

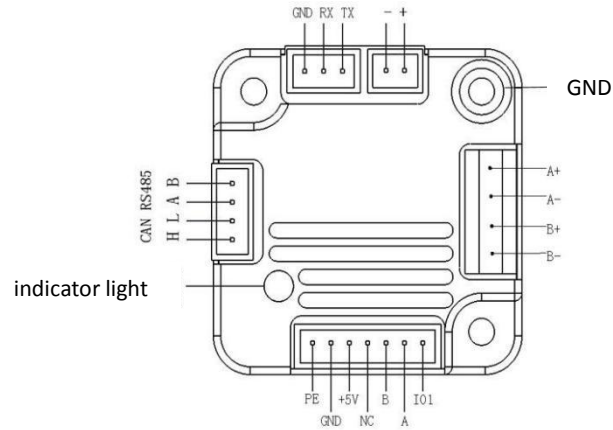

Port definition for driver control board:

Port	Description	Port	Description
+	DC24V Positive	B+, B-	Stepper motor Phase B wiring
-	DC24V Negative	IO ₁ (VM)	MOS positive
TX	RS232 Data input	IO ₂	IO2 signal
RX	RS232 Data output	IO ₃	IO3 signal
GND	RS232 Grounding	IO ₄	IO4 signal
A	RS485 A	+5V	Power positive
B	RS485 B	GND	GND
A+, A-	Stepper motor Phase A wiring	PE (OUT-)	MOS negative

Chapter 4 Software Communication

4.1 Brief Introduction

The data transmission between the syringe pump and the host controlling system (Computer, MCU, PLC, etc.) adopts serial communication (RS232, RS485). Description of the communication format as following: the communication adopts asynchronous serial communication, and the sum check with two bytes (2Byte) is adopted by the command & data frame. Commands & data in communication must be in hexadecimal. Parameters are stored in little-endian mode.

Communication interface: RS232, RS485;

Communication mode: two-way asynchronous, master-slave mode;

Baud rate: RS232&RS485: 9600bps, 19200bps, 38400bps, 57600bps, 115200bps;

Data bits: 8 Bit;

Parity check: no check;

Response time: <1 second after receiving the command

4.2 Installation & Debugging

1. Install the debugging tools, please refer to 《Debugging tools instructions》 for details.
2. Instructions for use, please refer to 《SY09S Quick Use Guide》 for details.

4.3 Communication Protocols

4.3.1 Command format

A: Interior parameters inquiry command

B: Device control command

C: Device interior parameter command setting

4.3.1.1 Common command format (Send 8 bytes, return 8 bytes)

Byte send:

B0	B1	B2	B3	B4	B5	B6	B7
FH (Frame header)	Address code	Function code	Status parameter		EOF (End of frame)	CUCUM (Cumulative sum)	
STX	ADDR	FUNC	1-8 Bit	9-16 Bit	ETX	Low byte	High byte

The 1st byte STX : Frame header (0xCC)

The 2nd byte ADDR : Slave address (0x00 ~ 0x7F)

Multicast address (0x80 ~ 0xFE) Broadcast address (0xFF)

- The 3rd byte FUNC : Function code
- The 4th, 5th byte : Parameters corresponding to the function code
- The 6th byte ETX : End of frame (0xDD)
- The 7th, 8th byte : Cumulative sum check code from byte 1 to 6

Byte return:

B0	B1	B2	B3	B4	B5	B6	B7
FH (Frame header)	Address code	Function code	Status parameter		EOF (End of frame)	CUCUM (Cumulative sum)	
STX	ADDR	STATUS	1-8 Bit	9-16 Bit	ETX	Low byte	High byte

- The 1st byte STX : Frame header (0xCC)
- The 2nd byte ADDR : Slave address (0x00 ~ 0x7F)
 Multicast address (0x80 ~ 0xFE) Broadcast address (0xFF)
- The 3rd byte FUNC : Function code
- The 4th, 5th byte : Parameters corresponding to the function code
- The 6th byte ETX : End of frame (0xDD)
- The 7th, 8th byte : Cumulative sum check code from byte 1 to 6

A : Interior parameters inquiry command (“Inquiry command”) (B2 ~ B4)

Command B2	Specification	Parameters B3、B4	Response Parameters B3、B4
0x20	Query address	0x0000	The value of address is 0x0000~0x007F and the default is 00.
0x21	Query RS232 baud rate	0x0000	Five baud rates: factory default is 9600bps. B3B4=0x0000 Corresponding baud rate is 9600bps. B3B4=0x0002 Corresponding baud rate is 38400bps. B3B4=0x0003 Corresponding baud rate is 57600bps. B3B4=0x0004 Corresponding baud rate is 115200bps.
0x22	Query RS485 baud rate		
0x27	Query maximum speed	0x0000	Maximum speed range of 3ml is 0x0000 ~ 0x0258. Maximum speed range of 8ml is 0x0000 ~ 0x01F4.
0x3F	Query current firmware version	0x0000	B3=0x01 B4=0x1E, as an example, indicate that the version is V1.30. See the version No. on the label for details.
0x4A	Query motor status	0x0000	Check status list for details
0x68	Query piston position	0x0000	B3B4=0x0000 After the pump is running, the command can be used to query the current position of the motor

			and display it The current distance between the motor and the zero position (in steps)
0x67	Synchronize piston position	0x0000	B3B4 = 0x0000 When the syringe pump is suddenly powered off during operation, the pump will continue to run for a short period of time, at which time the number of steps will be out of wrong, the number of steps to reach the reset position is not the home position when the power is turned on again. At this time, we need to run the position reset command 0x67, and then send 0x66 command to query current position, it means current position is the home position when B3=0x00 B4=0x00 received.
0x70	Query multicast channel 1 address	0x0000	B3B4 = 0x0000
0x71	Query multicast channel 2 address	0x0000	B3B4 = 0x0000
0x72	Query multicast channel 3 address	0x0000	B3B4 = 0x0000
0x73	Query multicast channel 4 address	0x0000	B3B4 = 0x0000

B : Device control command (“operation command”) (B2 ~ B4)

Command B2	Specification	Parameters B3、 B4	Response Parameters B3、 B4
0x42	Run in CW (clockwise)	Operating parameters	The value of B3&B4 ≥ 0 . When the number of steps corresponding to the B3&B4 parameter is greater than the maximum stroke steps, the motor will not run, and B3=02, B4=00 will return; when the number of steps is less than the Max. stroke steps, the devices will rotate according to the set steps
0x4D	Run in CCW (counterclockwise)	Operating parameters	The value of B3&B4 ≥ 0 . When the number of steps corresponding to the parameter B3&B4 is greater than the maximum number of stroke steps, the motor will not run and return B3=02, B4=00; when the number of steps corresponding to the parameter B3&B4 is less than the maximum number of stroke steps, the motor will follow the set number of steps.
0x4F (0x45)	Force reset	0x0000	B3B4 = 0x0000 When the syringe runs to the home position, the number of locked-rotor steps runs to the top through resetting which causes forced resetting. Then the piston will go back for a certain number of offset steps, leaving a little gap between

			the top of the piston and the syringe which greatly improve the service life of the piston seal. Note: command 0x4F forces a reset and does not recognize the dynamic speed set by 0x4B. 0x45 reset command, the speed of the first reset on power up defaults to 0x07 (the maximum speed set) and does not recognize the speed set by 0x4B (dynamic speed). To change the 0x45 reset speed, run 0x43 action once to recognize the dynamic speed set by 0x4B.
0x4B	Set dynamic speed (It will fail when power-off, and if not set, it will be the default speed.)	Operating parameters	The maximum speed of 3ml is 600rpm (0x0001 ~ 0x0258) . The maximum speed of 8ml is 300rpm (0x0001 ~ 0x012C) .
0x4E	Syringe to absolute position	Operating parameters	The absolute position of 3ml is 3600(0 ~ E10). The absolute position of 8ml is 3840(0 ~ F00).
0x49	Force stop	0x0000	B3B4 =0x0000
0x60	Open solenoid valve	B3B4 =0x0001	Switch to port NC
0x61	Close solenoid valve	B3B4 =0x0001	Switch to port NO

Status list

Status (B2)	Description of the response frame (B2) status
0x00	Normal status
0x01	Frame error
0x02	Parameter error
0x03	Optocoupler error
0x04	Motor busy
0x05	Motor stall
0x06	Unknown location
0x07	Command rejected
0x08	Illegal location
0xFE	Task pending
0xFF	Unknown error

Note: The code B2 in the response command indicates the current running status of the motor, only when B2=0x00, the motor runs normally, and the other parameters are as shown in the above table, which means different abnormal status respectively. In principle, the 0x4A command should be sent to query the motor status when the motor finishes operation. Only when the B2=0x00, other commands can be sent and execute correctly.

Caution: In RS485 communication, when sending an action command, the B2 byte in the response frame is FE, indicating that the command is received and is being executed.

4.3.1.2 Factory command format (send 14 bytes, return 8 bytes)

Factory commands need to be used with V0.8 debugging tools when RS232 or RS485 are used separately. See 《Quick Use Guide》 in detail.

Byte send:

B0	B1	B2	B3,B4 B5,B6	B7	B8	B9	B10	B11	B12	B13
FH (frame header)	Address code	Function code	Pass word	Function Parameter				EOF (End of frame)	CUCUM (Cumulative sum)	
STX	ADDR	FUNC	PWD	1-8 bit	9-16 bit	17-24 bit	25-32 bit	ETX	Low byte	High byte

The 1st byte STX : Frame header (0xCC)

The 2nd byte ADDR : Slave address (0x00 ~ 0x7F)

Multicast address (0x80 ~ 0xFE) Broadcast address (0xFF)

The 3rd byte FUNC : Function code

The 4th -7thbyte : Factory command format

The 8th -11thbyte : Parameters corresponding to the function code

The 12th byte ETX : End of frame (0xDD)

The 13th, 14thbyte : Cumulative sum check code from byte 1 to 12

Byte return:

B0	B1	B2	B3	B4	B5	B6	B7
FH (frame header)	Address code	Status code	Status Parameter		EOF (End of frame)	CUCUM (Cumulative sum)	
STX	ADDR	STATUS	1-8 Bit	9-16 Bit	ETX	Low byte	High byte

The 1st byte STX : Frame header (0xCC)

The 2nd byte ADDR : Slave address (0x00 ~ 0x7F)

Multicast address (0x80 ~ 0xFE) Broadcast address (0xFF)

The 3rd byte FUNC : Function code

The 4th-5thbyte : Parameters corresponding to the function code

The 6th byte ETX : End of frame (0xDD)

The 7th, 8thbyte : Cumulative sum check code from byte 1 to 6

C : Device interior parameter command setting ("Factory command")

Command B2	Abbreviation	Password B3 B4 B5 B6	Parameter Specification B7 B8 B9 B10
0x00	Set device address	B3=0xFF B4=0xEE B5=0xBB B6=0xAA	B7=0xFF (B8=0x00 B9=0x00 B10=0x00) The value range of XX is 00 ~ 7F and the default is 00.
0x01	Set RS232	B3=0xFF B4=0xEE	Totally 5 baud rates, the factory default is 9600bps.

	baud rate	B5=0xBB B6=0xAA	B7=0x0000 9600bps B7=0x0001 19200bps B7=0x0002 38400bps B7=0x0003 57600bps B7=0x0004 115200bps
0x02	Set RS485 baud rate	B3=0xFF B4=0xEE B5=0xBB B6=0xAA	
0x03	Set CAN baud rate	B3=0xFF B4=0xEE B5=0xBB B6=0xAA	Totally 4 baud rates, the factory default is 100K. B7=0x0000 corresponding baud rate 100Kbps B7=0x0001 corresponding baud rate 200Kbps B7=0x0002 corresponding baud rate 500Kbps B7=0x0003 corresponding baud rate 1Mbps
0x05	Set subdivision	B3=0xFF B4=0xEE B5=0xBB B6=0xAA	Totally 5 subdivisions: the factory default is subdivision 8. B7=0x0001 subdivision 2 B7=0x0002 subdivision 4 B7=0x0003 subdivision 8 B7=0x0004 subdivision 16 B7=0x0005 subdivision 32
0x07	Set maximum speed	B3=0xFF B4=0xEE B5=0xBB B6=0xAA	B7=0xFF (B8=0x00 B9=0x00 B10=0x00) The value range of FF is 0x01 ~ 0x258 and the default is 12C.
0x0E	Set automatic reset after power on	B3=0xFF B4=0xEE B5=0xBB B6=0xAA	B3B4=0x0000 do not automatically reset after power-on. B3B4=0x0001 automatically reset after power-on.
0x10	Set CAN destination address	B3=0xFF B4=0xEE B5=0xBB B6=0xAA	B7=0xFF (B8=0x00 B9=0x00 B10=0x00) The value range of FF is 00 ~ FF and the default is 00.
0x50	Set the multicast channel 1 address	B3=0xFF B4=0xEE B5=0xBB B6=0xAA	B7=0xFF (B8=0x00 B9=0x00 B10=0x00) The value range of FF is 80 ~ FE and the default value is 00.
0x51	Set the multicast channel 2 address	B3=0xFF B4=0xEE B5=0xBB B6=0xAA	B7=0xFF (B8=0x00 B9=0x00 B10=0x00) The value range of FF is 80 ~ FE and the default value is 00.
0x52	Set the multicast channel 3 address	B3=0xFF B4=0xEE B5=0xBB B6=0xAA	B7=0xFF (B8=0x00 B9=0x00 B10=0x00) The value range of FF is 80 ~ FE and the default value is 00.
0x53	Set the multicast channel 4 address	B3=0xFF B4=0xEE B5=0xBB B6=0xAA	B7=0xFF (B8=0x00 B9=0x00 B10=0x00) The value range of FF is 80 ~ FE and the default value is 00.

Example: Use the 0x50/51/52/53 command to set the multicast address (this example only uses the three commands 0x50/51/52)

Use three sets SY-09S syringe pumps with the same software version. In the RS485 communication mode, set their addresses to 00, 01, 02 and make a mark. Firstly, set parameter 0x81 of SY-09Smulticast channel 1 address marked as 00 to 81 by the command 0x50, set parameter 0x83 of multicast channel 3 address to 83 by command 0x52; Set parameter 0x81 of the multicast channel 1 address of SY-09S marked

as 01 to 81 by command 0x50, and the parameter 0x82 of multicast channel 2 address is set to 82 by command 0x51. Set parameter 0x82 of the multicast channel 2 address of SY-09S marked as 02 to 82 by command 0x51, and the parameter 0x83 of multicast channel 3 address to 83 by command 0x52 (See the table below for details).

Device Items	Device 1 (Address 0)	Device 2 (Address 1)	Device 3 (Address 2)
Multicast address	81	81	
		82	82
	83		83
Broadcast address	FF	FF	FF

After the setting is completed, connect the three devices in parallel to the serial port debugging tool, and use our debugging tool MotorTest V0.8 for debugging. In MotorTest V0.8, set the address to 0x81, the command to 0x4D and the parameter to 0xC8, click to send and then the pistons of device 1 and device 2 will have pumping action; Set the address to 0x82 ,the command to 0x4D and the parameter to 0xC8, click to send and then the pistons of device 2 and device 3 will have pumping action; Set the address to 0x83, the command to 0x4D and the parameter to 0xC8, click send and then the pistons of device 1 and device 3 will have pumping action; Set the address to 0xFF, the command to 0x4D, and the parameter to 0xC8 , click to send and then the pistons 1, 2, and 3 will all have pumping action.

The newly added command to set the multicast address meets the needs of customer groups to a great extent and makes it easier for customers to select the equipment that you want to control, and can complete the work requirements more efficiently and quickly during use.

Note:

All the parameters of all the above commands are set in little-endian mode. In little-endian mode storage, the low bit of data is stored in the low bit of the address, and the high bit of data is stored in the high bit of the address.

Chapter 5 Common Problems & Solutions

Item	Fault	Reason	Troubleshooting method
1	Not working	The working voltage is not in the acceptable range.	Check whether the actual voltage deviates from the rated voltage.

	when powered on	The connection is loose or disconnected.	Manually check whether the connection is good, or check the line with a multimeter.
2	Unable to aspirate or aspirate properly	The pipe system is not tightly sealed.	Check whether the joint is tight.
		The aspirating pipe is blocked.	Clean and dredge the pipe.
		The aspirating valve or the dispensing valve is blocked by debris.	Clear the debris.
3	Bubbles	Air leakage in aspirating pipe.	Find the leak and eliminate it.
		The inlet and outlet pipe joints are not tightly sealed.	Replace the gasket and tighten the pipe joint.
		Gasket broken	Replace the gasket.
		Excessive fluid pipe diameter variation	The diameter of the fluid path should be as consistent as possible.
4	Pump stuck	Optocoupler is not triggered.	Check the optocoupler wiring (refer to the optocoupler wiring method).
		Optocoupler is burned out.	Replace optocoupler.
		Reverse connection of motor wires.	Switch any phase of the motor wires.
5	Motor overheated	Drive voltage is too large.	Adjust voltage
		Drive current is too large.	Adjust current
		Holding current is too large.	Holding current $\leq 50\%$ of rated current
6	Abnormal sound of pump operation	Motor running speed is too high or too low.	Adjust the motor speed to a suitable value.
		There are crystals in the pump head.	Cleaning steps are required after the machine runs or before it starts.
7	Poor sampling accuracy	The pipe system is not tightly sealed.	Check whether the joint is tight.
		There are bubbles in the pipe.	Refer to troubleshooting method of above point 2.

Chapter 6 Quick Command

A Quick Commands

Code B2	Abbreviation	Parameter Description B3 B4
0x20	Query address	Address ranging from: 0x0000 ~ 0x007F, default 00
0x27	Query the maximum speed	5ml, 12.5ml maximum speed range is 0x0000 to 0x0258 the maximum speed for 25ml is 0x0000 to 0x01F
0x3F	Query current version	B3=0x01 B4=0x09, above is just an example, if the response parameter is the same as above parameter, it means the current version is V1.9, see the version number on the label for details

0x4A	Query motor status	B3=0x00 B4=0x00
0x4D	Aspirate liquid	When the number of steps corresponding to parameter B3B4 is greater than the maximum number of steps, the motor will not run, and return byte B3=08, B4=00; When the number of steps corresponding to B3B4 parameter is set to be less than the maximum number of steps, the motor rotates according to the set number of steps.
0x42	Discharge liquid	When the number of steps corresponding to parameter B3B4 is greater than the maximum number of steps, the motor will not run, and return byte B3=08, B4=00; When the number of steps corresponding to B3B4 parameter is set to be less than the maximum number of steps, the motor rotates according to the set number of steps.
0x45	Reset of syringe pump	B3=0x00 B4=0x00 The syringe pump runs to the home position and stops
0x4F	Forced reset of syringe pump	B3=0x00 B4=0x00 The switching valve runs to the encoder home position, which overlaps with the position reset by the 0x45 command
0x49	Force Stop	B3=0x00 B4=0x00
0x4B	Set dynamic speed	B3B4 ranges from 0x0001 to 0x01C2. The speed of motor is 1 ~ 450
0x4E	Syringe runs to absolute position	B3B4 ranges from 0x0000 ~ 0x1770, which is optional position in the syringe stroke
0x60	Valve open	Switch port to NC port Note: When the environment and fluid temperature of rubber membrane materials FKM and Kalrez are below 15 C° The switching time of the valve is extremely long, please pay attention to it.
0x61	Valve closed	Switch port to NO port (default) Note: When the environment and fluid temperature of rubber membrane materials FKM and Kalrez are below 15 C° The switching time of the valve is extremely long, please pay attention to it.

B Hexadecimal Quick Reference Table

Decimal	Hexadecimal	Decimal	Hexadecimal	Decimal	Hexadecimal	Decimal	Hexadecimal
1	1	110	6E	300	12C	1000	3E8
5	5	120	78	325	145	1100	44C
10	A	130	82	350	15E	1200	4B0
15	F	140	8C	375	177	1300	514
20	14	150	96	400	190	1400	578
25	19	160	A0	425	1A9	1500	5DC
30	1E	170	AA	450	1C2	1600	640
35	23	180	B4	475	1DB	1700	6A4
40	28	190	BE	500	1F4	1800	708
45	2D	200	C8	525	20D	1900	76C

50	32	210	D2	550	226	2000	7D0
55	37	220	DC	575	23F	2100	834
60	3C	230	E6	600	258	2200	898
65	41	240	F0	650	28A	2300	8FC
70	46	250	FA	700	2BC	2400	960
75	4B	260	104	750	2EE	2500	9C4
80	50	270	10E	800	320	2600	A28
90	5A	280	118	850	352	2700	A8C
100	64	290	122	900	384	2800	AF0

C Switching Protocol

1. SY08 has two protocols, RUNZE and ASCII protocol. If you need to switch protocols, you can refer to the following steps (both need to be powered off after switching)

- Burn-in Protocol
- If the default programmed command is the RUNZE command, it will query the following return code

Send→ 91 EB 07 00 00 00 00 00 D5 28 FF F8

Receive← 91 EB 02 01 00 63 D7 F6 AB 00

2. parameter definition: 02 stands for RUZNE

- If the default programmed command is the ASC II command, it will query the following return code

Send→ 91 EB 07 00 00 00 00 00 D5 28 FF F8

Receive← 91 EB 0A 01 00 02 C4 47 0B 00

0A parameter definition: 0A stands for ASC II

- Switching protocol (must use RS232 port to switch the address, and must be powered off after switching)
- If RUNZE command is required, the following code can be used to switch the address

Set the address to 0x0802 (RUNZE)

Send→ 91 EB 03 00 00 02 08 00 00 0C 0A 69 69

Receive← 91 EB 00 01 00 0D 5A 8A 40 00

The values queried after the setting are as follows.

Send→ 91 EB 07 00 00 00 00 00 D5 28 FF F8

Receive← 91 EB 02 01 00 63 D7 F6 AB 00

- If ASC II command is required, the following code can be used to switch the address

Send→ 91 EB 03 00 00 0A 08 00 00 6D 19 D8 C9

Receive← 91 EB 00 01 00 0D 5A 8A 40 00

The values queried after the setting are as follows.

Send→ 91 EB 07 00 00 00 00 00 00 D5 28 FF F8

Receive← 91 EB 0A 01 00 02 C4 47 0B 00

Chapter 7 Version Description

Version	Description	Release Time
V1.0	Initial version	2023.3.28
V1.1	Add precautions for solenoid valves	2024.1.30
V1.2	Modify the command to query the syringe pump	2024.4.2

Chapter 8 Technical Service

Nanjing Runze Fluid Control Equipment Co.,LTD

Landline (FAX)	025-5119 7362
SALE	+86 173 6638 4502
Technical Service	+86 198 2581 4316
Email	runzeliuti@runzeliuti.com
Website	www.runzeliuti.com
Shop	https://runzeliuti.en.alibaba.com
Address	No.9 Tianxing West Road, Dongshan Street, Jiangning District, Nanjing, Jiangsu Province, China



Official URL



Alibaba Store URL



Aliexpress Store URL